

GelNest™ Matrix, Growth Factor Reduced

Gel Concentration Inquiry: If the label on the bottle is lost, please click or copy the link below to your browser, and select "Related Reference Tools-> Certificates" to download the batch-specific COA.

<https://www.nestscientificusa.com/product/detail/636882319776419840>

Product overview

GelNest™ Matrix is prepared from basement membrane components extracted from mouse tumor tissues. The main components include laminin, type IV collagen, heparan sulfate proteoglycan, etc. These components can provide the support and signals required for cell adhesion, differentiation, and proliferation. They can also simulate the characteristics of the basement membrane in a physiological environment and improve the success rate and effect of cell culture.

In addition to basement membrane components, GelNest™ Matrix is also rich in a variety of growth factors. These growth factors can promote cell differentiation, proliferation, and migration, further mimicking cell signaling pathways and interactions in physiological environments. GelNest™ Matrix has a wide range of application prospects, especially in tissue engineering, cell culture and research. It can be used for research on organoid culture, stem cell differentiation, angiogenesis, migration or invasion, and *in vivo* tumorigenesis.

Product information

Product number	Product name	Packaging specifications
211232	GelNest™ Matrix, Growth Factor Reduced, LDEV-Free	Bag Package, 5 mL/bottle, 1 bottle/bag
211242	GelNest™ Matrix, Growth Factor Reduced, Phenol Red-Free, LDEV-Free	Bag Package, 5 mL/bottle, 1 bottle/bag

211231	GelNest™ Matrix, Growth Factor Reduced, LDEV-Free	Bag Package, 5 mL/bottle, 2 bottle/bag
211241	GelNest™ Matrix, Growth Factor Reduced, Phenol Red-Free, LDEV-Free	Bag Package, 5 mL/bottle, 2 bottle/bag
211332	GelNest™ Matrix, Growth Factor Reduced, LDEV-Free, Ultra-low Endotoxin	Bag Package, 5 mL/bottle, 1 bottle/bag
211342	GelNest™ Matrix, Growth Factor Reduced, Phenol Red-Free, LDEV-Free, Ultra-low Endotoxin	Bag Package, 5 mL/bottle, 1 bottle/bag

Product parameters

Source	Mouse tumor tissue basement membrane components
Formulation*	Growth factor reduced
Protein concentration	See label, or please download the COA from our official website to obtain a lot-specific concentration.
Appearance	GelNest™ Matrix is liquid at 4°C but forms a gel at 37°C. Phenol red-containing gel appears bright yellow when frozen, and red at temperatures above 0° C.
Applications	Suitable for 2D primary cell culture, organoid construction, culture, differentiation, invasion assays
Storage and shelf life	Store in a refrigerator at -20°C (frost-free function off) or a -80°C freezer for up to 2 years. It is recommended to aliquot the thawed product into single-use portions and store it in -20°C or -80°C for up to 2 years.
Precautions	GelNest™ Matrix will start to solidify when the temperature is higher than 10°C. Please try to operate on ice as much as possible, and it is recommended to pre-cool the consumables that directly contact the gel, such as pipette tips.

*Please use phenol red-free matrix gel for colorimetric analysis.

Experimental procedures

GelNest™ Matrix Coating Recommendations

Application Type	Dilution Factor	Final Concentration	Volume Added	Experiments
Thin Gel	1:10	>1mg/mL	50µL/cm ²	2D primary cell culture
	1:50	>0.1mg/mL		Invasion assay with cell culture inserts
	1:100	>0.01mg/mL	300µL/cm ²	Stem cell culture
Thick Gel	Gel : Cell Mixture<7:3	>7mg/mL	150-200µL/cm ²	Organoid culture
	No Dilution	>10mg/mL		Angiogenesis

Please determine the specific experimental steps based on cell types, culture conditions, and application experience.

Organoid culture

1. Re-suspend the single cell suspension used for organoid culture in pre-cooled basal medium at 4°C, and count the cells.
2. Mix GelNest™ Matrix with the cells (dilution factor<70% recommended) and add the mixture to a preheated 24-well plate, each well containing approximately 5x10⁴ cells and 60µL of the matrix gel.
3. Immediately place the well plate into the incubator. After about 10 minutes, the matrix gel will solidify.
4. Add 500µL of organoid culture medium for culture.
5. Wait 3-5 days for the organoids to form. Finally, the sensitivity of organoids to various drugs can be determined by imaging live cells through high-content microscopy.

* It is recommended to use GelNest™ Matrix, for Organoid Culture (NEST 211282) for better results.

Invasion experiment

1. Use HT-1080 cells in MEM medium supplemented with 10% fetal bovine serum and culture them to a cell confluence of 80% to 90% before use.
2. Take 20 μ L of GelNest™ Matrix, dilute it to 1000 μ L with serum-free MEM (1:50 dilution).
3. Gently pipette up and down to thoroughly mix the gel.
4. Add 100 μ L of the diluted matrix gel mixture to the center of each cell culture insert (NEST transwell product) so that the matrix gel mixture evenly covers the surface of the insert. Incubate the culture dish at 37°C for 1 hour to allow gel formation.
5. After trypsinization of the cells (typically, for a 6-well plate, digest the cells with 200 μ L of trypsin at 37°C for 3 minutes, then terminate the digestion with 10% serum, centrifuge at 300xg for 3 minutes), resuspend the cells in serum-free MEM culture medium.
6. Count cells and take 750 μ L of cells at a starting density of 1x10⁶/mL (expected to use 10 wells with 7.5x10⁴ cells per well, a total of 750,000 cells), and dilute it with MEM serum-free medium to 1.5mL.
7. Seed 150 μ L of cell suspension into the upper chamber of each cell culture insert, resulting in 7.5x10⁴ cells/well.
8. In the experimental group, add 800 μ L of culture medium containing 10% FBS as a chemoattractant to the lower chamber, while in the control group, add 800 μ L of serum-free culture medium to the lower chamber. Incubate the cells overnight at 37°C with 5% carbon dioxide in a humidified incubator.
9. Discard the supernatant medium from the cell culture insert and wash twice with PBS. Stain the cells on the underside of the membrane with crystal violet for 10 minutes.
10. Wash the cell culture inserts twice with PBS to remove unbound crystal violet. Use a moist cotton swab to gently remove the cells from inside the cell culture insert, then air-dry.
11. Observe and capture images of the invaded cells under a microscope.
12. Dilute acetic acid to 33% (v/v) with ddH₂O to elute the bound crystal violet. Add

400µL 33% acetic acid to each cell culture insert and shake on a shaker for 10 minutes.

Transfer the eluate from the lower chamber to a 96-well transparent microplate, and

measure the absorbance at 590 nm using a microplate reader.

Safety recommendations and limitations

Please follow good laboratory safety practices.

For research use only. Not intended for diagnostic or therapeutic purposes. Contains ingredients of animal origin.

Technical support and contact information

For FAQ, GelNest™ Matrix Selection Guide, Quality Assurance COA/COC or other technical support and product issues, please refer to our website or use the following contact information.

Production and after-sales service unit: Wuxi NEST Biotechnology Co., Ltd.

Production and after-sales address: No.530, Xida Road, Meicun Industrial Park, Xinwu District, Wuxi, Jiangsu, China

Tel: +86-510-68006788

Email: info@nest-wuxi.com

Website: www.cell-nest.com